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## Photobiological Testing Instrument

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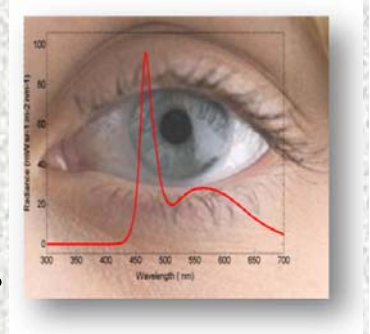
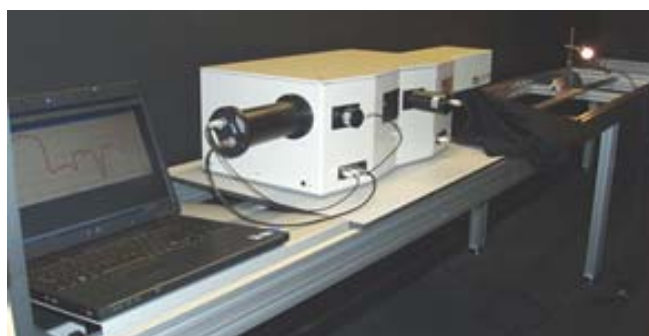
Apparent source location:		35mm behind plane of front face																
Apparent size of the array:		177 mrad																
Apparent size of component LED:		15x15 mrad																
<b>EN62471:2008 Evaluation</b>																		
Measured at 200mm from apparent source																		
Irradiance-based hazards																		
Hazard	Measured value	Allotted risk group	Risk group limit value	Time to exposure limit (s)														
Actinic UV ( $mW m^{-2}$ )	6.0	RG2	30	5025														
Near UVA ( $W m^{-2}$ )	62.8	RG2	100	159														
Radiance-based hazards																		
Hazard	Measured value ( $W sr^{-1} m^{-2}$ )	Risk group limit value ( $W sr^{-1} m^{-2}$ )	Risk group tested	Pass/Fail														
Blue light hazard 100mrad FOV	37.1	100	Exempt	Pass														
Retinal Thermal 11mrad FOV	$3.94 \times 10^6$	$1.87 \times 10^6$	Exempt	Pass														
Spectral Irradiance at 200mm		Overview of EN62471:2008 classification																
		<table border="1"> <thead> <tr> <th>Hazard</th> <th>Risk Group</th> </tr> </thead> <tbody> <tr> <td>Actinic UV</td> <td>RG2</td> </tr> <tr> <td>Near UVA</td> <td>RG2</td> </tr> <tr> <td>Blue Light</td> <td>Exempt</td> </tr> <tr> <td>Retinal Thermal</td> <td>Exempt</td> </tr> <tr> <td>IR eye</td> <td>Exempt</td> </tr> <tr> <td>Thermal Skin</td> <td>Exempt</td> </tr> </tbody> </table>			Hazard	Risk Group	Actinic UV	RG2	Near UVA	RG2	Blue Light	Exempt	Retinal Thermal	Exempt	IR eye	Exempt	Thermal Skin	Exempt
Hazard	Risk Group																	
Actinic UV	RG2																	
Near UVA	RG2																	
Blue Light	Exempt																	
Retinal Thermal	Exempt																	
IR eye	Exempt																	
Thermal Skin	Exempt																	
IEC62471-2 Labelling																		
Risk group 2 Caution: UV emitted from this product. Minimise exposure to eyes or skin. Use appropriate shielding.																		
AORD Evaluation																		
Measured at point of closest proximity																		
Hazard	Measured Value	Allotted Risk Group	Risk Group Limit Value	Time to Exposure Limit (s)														
Actinic UV ( $mW m^{-2}$ )	65.3	RG3	-	480														
Near UVA ( $W m^{-2}$ )	686.1	RG3	-	15														

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Bentham Instruments Limited